
DCAT 15
The dynamic contact angle measuring instrument and tensiometer



**dataphysics** 

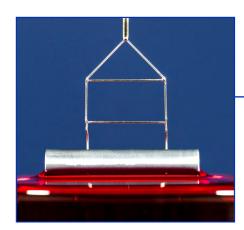

Density determination of liquids with the density determination set DIS 11


The dynamic contact angle measuring instrument and tensiometer **DCAT 15** is the standard instrument for the weight-based measurement of contact angle, surface and interfacial tension, critical micelle formation concentration, density, sedimentation and penetration.



holders for fibre bundles FH 11 and powders PUR 11



film holder FO 11 and plate holder PSH 11

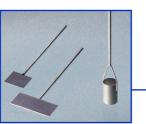

Wilhelmy plates PT 11/9 and cylindrical plate PT 10



penetration probe PP 11 and sedimentation cone SC 11

Software for efficient work

for the DCAT 15 are:

The newly developed, Windows® based,

DCATS 31 — surface/interfacial tension

• determination of the surface and inter-

DCATS 32 — dynamic contact angle

• determination of the dynamic contact

angle of solids (e.g. plates, films, rods

• adsorption measurement on powders

 analysis of the surface free energy of solids as well as its components accord-

ing to nine different theories


facial tension according to the Du Noüy ring and the Wilhelmy plate method

discretely usable modules, and is operable traditionally, using mouse and keyboard,

**DCATS**oftware is available in various




density determination set for solids DSS 11/12

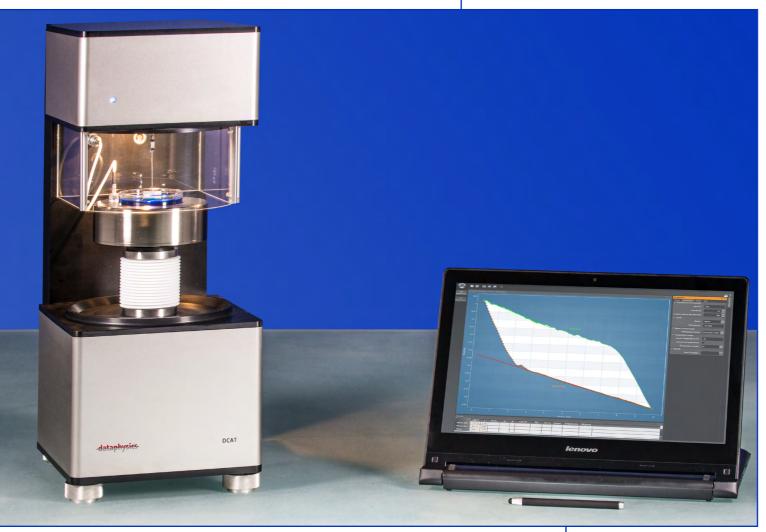

Determination of the contact angle hysteresis with a DCAT 15 with Wilhelmy plate PT 11

## Main features of the DCAT 15

- high-precision electrodynamic compensation weighing system with automatic and manual calibration
- software controlled, motor-driven height positioning of the sample receptacles with variable speed
- $\bullet$  automatic coupling lock for the balance
- automatic crash protection for measuring probes during measurement
- illuminated sample chamber
- integrated magnetic stirrer
- digital thermometer with connections for two Pt 100 temperature sensors

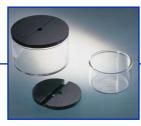

sample vessels GS 70/50 and cover plates CP 50/70



density determination set DIS 11 and Du Noüy rings RG 11/10

# or on multi-touch notebooks/pads by finger/pen. The available software modules

• determination of the density of liquids

## DCATS 35 — sedimentation/penetration

 determination of the sedimentation rate, yield forces, the penetration resistance and rate

## DCATS 36 — solid density

DCATS 34 — liquid density


• determination of the density of solids

### DCATS 33 — CMC

and fibres)

and fibre bundles

 automated determination of the critical micelle formation concentration (CMC) of surfactants, using the dosing unit LDU x/x for additive and subtractive dosing

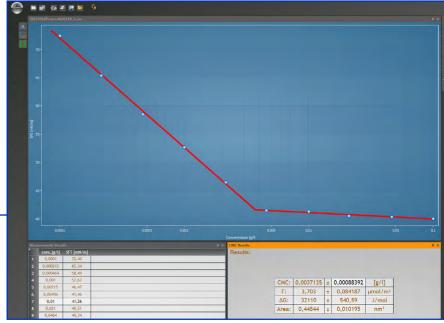

DCATS 33 — determination of the critical micelle formation concentration



LDU 2/2 dosing system for the automated CMC determination

# fechnical information in this document is subject to change without prior notice. Errors and omissions excepted. © Copyright by DataPhysics Instruments GmbH, Filderstadt. dataphysics is a registered trademark of DataPhysics Instruments GmbH DS/DCAT15-1604 – 1.0/Fin Photos by Norbert Heil, Thomas Müller, Bavaria, Tony Stone, Dreamstime, DataPhysics archives. Artwork and layout: Daniel Maier

# **dataphysics**

# Technical data

| Measuring range for surface and interfacial tensions: | • 1 1000 mN/m; ± 0.01 mN/m resolution  |
|---|--|
| Measuring range for contact angles:                   | • 0 180°; ± 0.01° resolution   |
| Measuring range for densities:                        | • 0.50 2.50 g/cm³; ± 0.002 g/cm³ resolution  |
| Weighing range:                                       | • 100 µg 220 g   |
| Measuring value range:                                | • up to 50 weighing values per second  |
| Traversing range and speed for sample table:          | • 80 mm<br>• 46 nm/s 12 mm/s   |
| Travel resolution:                                    | • 24 nm  |
| Balance calibration:                                  | • automatic internal and manual external with reference weights  |
| Automatic stirrer:                                    | • integrated, software controlled  |
| Temperature measurement and range:                    | <ul> <li>-10 130 °C (liquid temperature control unit TV xx)</li> <li>2 x Pt 100 inputs for -60 +450 °C (Pt 100 as option); ± 0.01 K resolution; precision 1/3 DIN IEC 751 (±0.03%), Class B</li> </ul> |
| Dimensions (L x W x H):                               | • 250 x 205 x 500 mm <sup>3</sup>  |
| Weight:   | • 15 kg  |
| Power supply:   | • 100 240 VAC; 50 60 Hz; 70 W  |

# Standards

The high degree of accuracy of the DCAT devices complies with all related international standards, for example:

- $\bullet \ \textbf{ISO 6295} \ \text{Petroleum products} \text{Mineral oils} \text{Determination of interfacial tension of oil against water} \text{Ring method}$
- ISO 6889 Surface active agents; Determination of interfacial tension by drawing up liquid films
- ASTM D971 Standard Test Method for Interfacial Tension of Oil Against Water by the Ring Method
- **ASTM D1417** Standard Test Methods for Rubber Latices-Synthetic
- DIN EN 14210 Surface active agents Determination of interfacial tension of solutions of surface active agents by the stirrup or ring method
- ASTM D1331 Standard Test Methods for Surface and Interfacial Tension of Solutions of Paints, Solvents, Solutions of Surface-Active Agents, and Related Materials
- ISO 304 Surface active agents; Determination of surface tension by drawing up liquid films
- DIN ISO 1409 Plastics/rubber Polymer dispersions and rubber latices (natural and synthetic) Determination of surface tension by the ring method
- $\bullet \ \ \textbf{OECD 115} \ \ \textbf{OECD Guidelines for the Testing of Chemicals: Surface Tension of Aqueous Solutions}$

# Accessories (excerpt)

set of reference weights **RWS** • set of reference weights, DKD certified **RWS-C** • liquid temperature control unit for sample vessels with diameters of 50 mm (**TV 50**) and 70 mm (**TV 70**) with integrated Pt100 probe • automatic dosing and refill system **LDU xx** • sample vessels made of glass **GS xx** and PTFE **GS xxP** as well as cover plates **CP xx** • Du Noüy rings **RG 11/10** • Aligning tool **R-AT** • Wilhelmy plates **PT 11/9** • cylindrical plate **PT 10** • density determination set for liquids **DIS 11** • density determination set for solids **DSS 11/12** • sedimentation cone **SC 11** • penetration probe **PP 11** • kit for measurement of powder samples **PUR 11** • holder for powders, pigments, fibres and fibre bundles **FH 11** • plate holder **PSH 11** • film holder **FO 11** • filter papers for PUR 11 **FP 11** and FH 11 **FP 12** 

For more information about a tailor-made solution to your surface chemistry requirements, please contact us.

We will be pleased to provide a quotation, obligation free, for your instrument system.

Your sales partner:

DataPhysics Instruments GmbH • Raiffeisenstraße 34 • 70794 Filderstadt, Germany phone +49 (0)711 770556-0 • fax +49 (0)711 770556-99 sales@dataphysics.de • www.dataphysics.de